


ProScope 5Mp Microscope Camera - QuickStart Guide

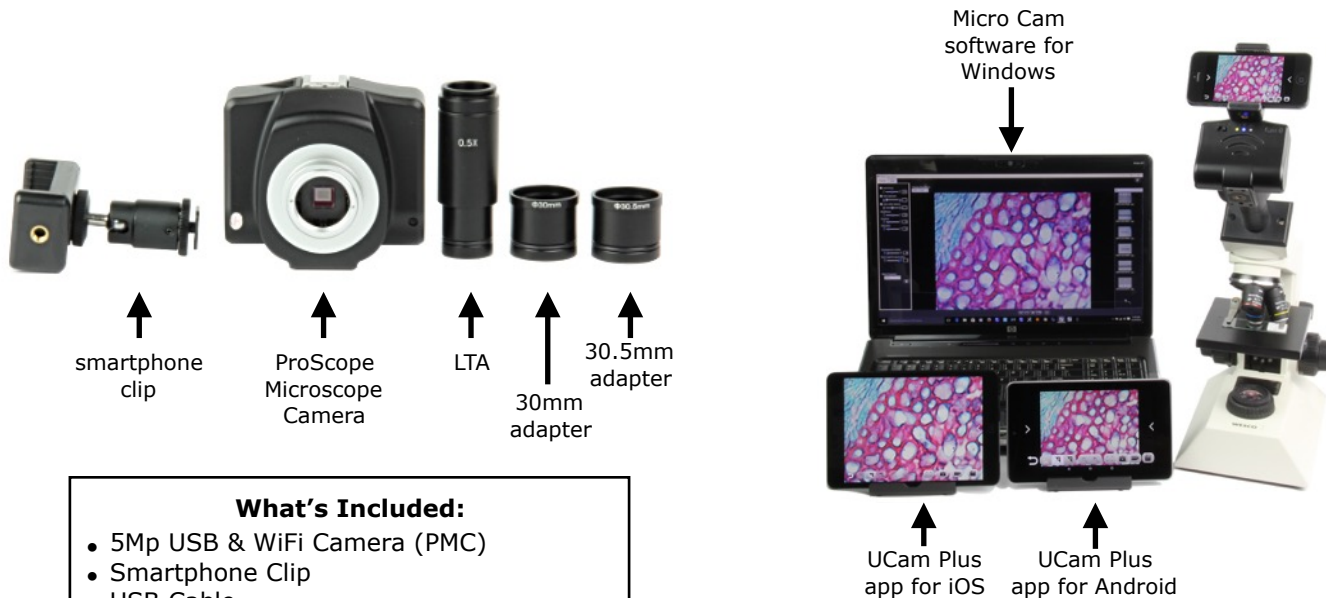

What's Included:

- 5Mp USB & WiFi Camera (PMC)
- Smartphone Clip
- USB Cable
- USB Charging Adaptor
- 23mm Lens Tube Adaptor (LTA)
- 30mm microscope tube adapter
- 30.5mm microscope tube adapter
- Windows Software on CD
- iOS app - [UCam Plus](#) free download
- Android app - [UCam Plus](#) free download
- Mac Software - free download of [ProScope HR software](#)
- Chrome OS Software - use native camera app or [Webcam Toy](#)

SPECIFICATIONS

Model #: PS-MC5UW
Description: 5Mp USB & WiFi Digital Eyepiece Camera for microscopes
Imager: 5Mp Aptina color CMOS
Interface: USB & WiFi
Desktop Software Included: Windows
Apps included: iOS, Android
WiFi compatibility: Windows, iOS, Android
USB compatibility: Windows, Mac, Chrome
Regulatory compliance: CE, FCC, RoHS
Simultaneous connected WiFi devices: 7 max
Power Source: internal battery or 5v USB
Battery life: approximately 8 hrs

Getting Started

1. Charge the internal battery of the ProScope Microscope Camera (PMC) with the included USB cable and power adapter, or from the USB port of your computer.
2. For WiFi operation, turn on the PMC by selecting the WiFi switch setting on the back of the unit.
3. Attach the LTA to the PMC and slide the LTA into your microscope's receiving tube after removing the microscope's eyepiece. Use the provided 30mm or 30.5mm adapters if needed.
4. For Windows:
Install the Micro Cam software from the included CD.
From your network control panel, choose the VideoWiFi-40xxxx network and enter the default password 12345678.
Launch the Micro Cam software and accept using the WiFi camera as prompted.
5. For iOS (iPad, iPhone):
Download the free UCam Plus app from the iTunes App Store.
From your device's Settings, choose the VideoWiFi-40xxxx network and enter the default password 12345678.
Launch the UCam Plus app and tap on the preview icon to view a live image from the PMC.
6. For Android OS:
Download the free UCam Plus app from the Google Play App Store.
From your device's Settings, choose the VideoWiFi-40xxxx network and enter the default password 12345678.
Launch the UCam Plus app and tap on the preview icon to view a live image from the PMC.
7. For USB operation, connect the provided USB cable to the PMC then to the USB port on your computer.
Turn on the PMC by selecting the USB switch setting on the back of the unit. Note: USB and WiFi modes do not operate simultaneously. For Windows, launch the Micro Cam software and select USB camera operation. For Mac and Chrome operation, see options listed in the What's Included section above.