

ProScope HR Quick Start

Macintosh:

1. Make sure you have Mac OS 10.4.8 and QuickTime 7 or later installed on your computer.*
2. Launch (double-click) the ProScope HR software installer located on the included CD-ROM. Macintosh folder>ProScope HR Installer.
3. Please quit iChat and unplug (or turn off) any external video camera before plugging in the ProScope HR USB microscope.
4. Plug the ProScope HR USB microscope into a USB 2.0 high speed port on your computer.**
5. Launch (double-click) the ProScope HR software application located in your applications folder. Now you can begin using your ProScope HR microscope.***

Windows:

1. Make sure you have Windows XP Service Pack 2 (also Windows Vista compatible) and QuickTime 7 or later installed on your computer.*
2. Launch (double-click) the ProScope HR software installer located on the included CD-ROM. Windows folder>InstallProScopeHR.
3. Plug the ProScope HR USB microscope into a USB 2.0 high speed port on your computer.**
4. Launch (double-click) the ProScope HR software application from your desktop ProScope HR icon or "Programs" Bodelin folder. Now you can begin using your ProScope HR microscope.***

**QuickTime is included on the CD-ROM for your convenience. The Windows ProScope HR installer will automatically install QuickTime on your computer if necessary.*

***Note – there is no driver installation required since the ProScope HR microscope uses the native OS USB Video Class and HID Class driver available in Mac OS 10.4 and Windows XP Service Pack 2 or later.*

****Make sure you plug in the ProScope HR USB microscope BEFORE launching the ProScope HR software application.*

Important READ ME FIRST files and additional documentation are available on the CD-ROM provided or online at www.proscopehr.com.